

Chromium Connect Software Release Note

FOR USE WITH

Chromium Connect, PN-1000171

Chromium Connect, PN-1000180

Notices

Document Number

CG000352 • Rev C

Legal Notices

© 2020 10X Genomics, Inc (10x Genomics). All rights reserved. Duplication and/or reproduction of all or any portion of this document without the express written consent of 10x Genomics, is strictly forbidden. Nothing contained herein shall constitute any warranty, express or implied, as to the performance of any products described herein. Any and all warranties applicable to any products are set forth in the applicable terms and conditions of sale accompanying the purchase of such product. 10x Genomics provides no warranty and hereby disclaims any and all warranties as to the use of any third-party products or protocols described herein. The use of products described herein is subject to certain restrictions as set forth in the applicable terms and conditions of sale accompanying the purchase of such product. A non-exhaustive list of 10x Genomics' marks, many of which are registered in the United States and other countries can be viewed at: www.10xgenomics.com/trademarks. 10x Genomics may refer to the products or services offered by other companies by their brand name or company name solely for clarity, and does not claim any rights in those third party marks or names. 10x Genomics products may be covered by one or more of the patents as indicated at: www.10xgenomics.com/patents. All products and services described herein are intended FOR RESEARCH USE ONLY and NOT FOR USE IN DIAGNOSTIC PROCEDURES.

Instrument & Licensed Software Updates Warranties

Updates to existing Instruments and Licensed Software may be required to enable customers to use new or existing products. In the event of an Instrument failure resulting from an update, such failed Instrument will be replaced or repaired in accordance with the 10x Genomics Limited Warranty, Assurance Plan or service agreement, only if such Instrument is covered by any of the foregoing at the time of such failure. Instruments not covered under a current 10x Genomics Limited Warranty, Assurance Plan or service agreement will not be replaced or repaired.

Support

Email: support@10xgenomics.com

10x Genomics

6230 Stoneridge Mall Road

Pleasanton, CA 94588 USA

Chromium Connect Software Release Note

Item	Description
For Use With	PN-1000171 Chromium Connect PN-1000180 Chromium Connect
Software Version	V2020.10.07
Software Release Date	October 2020
Compatibility	Chromium Connect Instrument
Features	<ol style="list-style-type: none">1. Loading of pipette tip racks scales with number of samples2. LIMS Output to include a run result field (i.e. SUCCESS FAILURE ABORT)3. Screen display to indicate status of Chromium Connect Software update
Bug Fixes & Changes	<ol style="list-style-type: none">1. HEPA light remains ON during a run2. cDNA no longer aliquoted to TS-1 when mid-run QC is skipped3. Updated maintenance screens4. Pre-run deck check ensures Waste Bar is present5. Repetitive prompt to refill pipette tips removed6. Wraparound selection of SI wells now compatible with the assay7. Secondary Assay transition corrected to proceed with beginning or end of run inputs

Previous Software Releases

Previous Software Version V2020.07.31
Released: August 2020

Features

1. Chromium Automated Controller Firmware update (v4.1.0) integrated into Chromium Connect Software update
2. Software integration for HEPA CAP hood functionality
3. Capability to export barcode information into LIMS and run audit logs
4. Added LIMS compatibility for qPCR and Pooling

Bug Fixes & Changes

1. Updated User Interface screens for general consistency of language and formatting
 2. Added sound notification for GEM Check Timer
 3. Minor bug fixes to user interface and barcode scanning
 4. Added ability to remove network drives in menu settings
 5. Enabled data purging off storage databases
 6. Easier access to waste bin during maintenance
 7. Pooling assay accepts sample input
-